


# MGMR<sup>®</sup> Versatile Nano-vehicle for Non-viral Cellular Transfection

Aaron Tasset, Kevin Castillo, Milos Marinkovic


Biopact, LLC

12310 Trail Driver St. 78737

Austin, Texas, USA

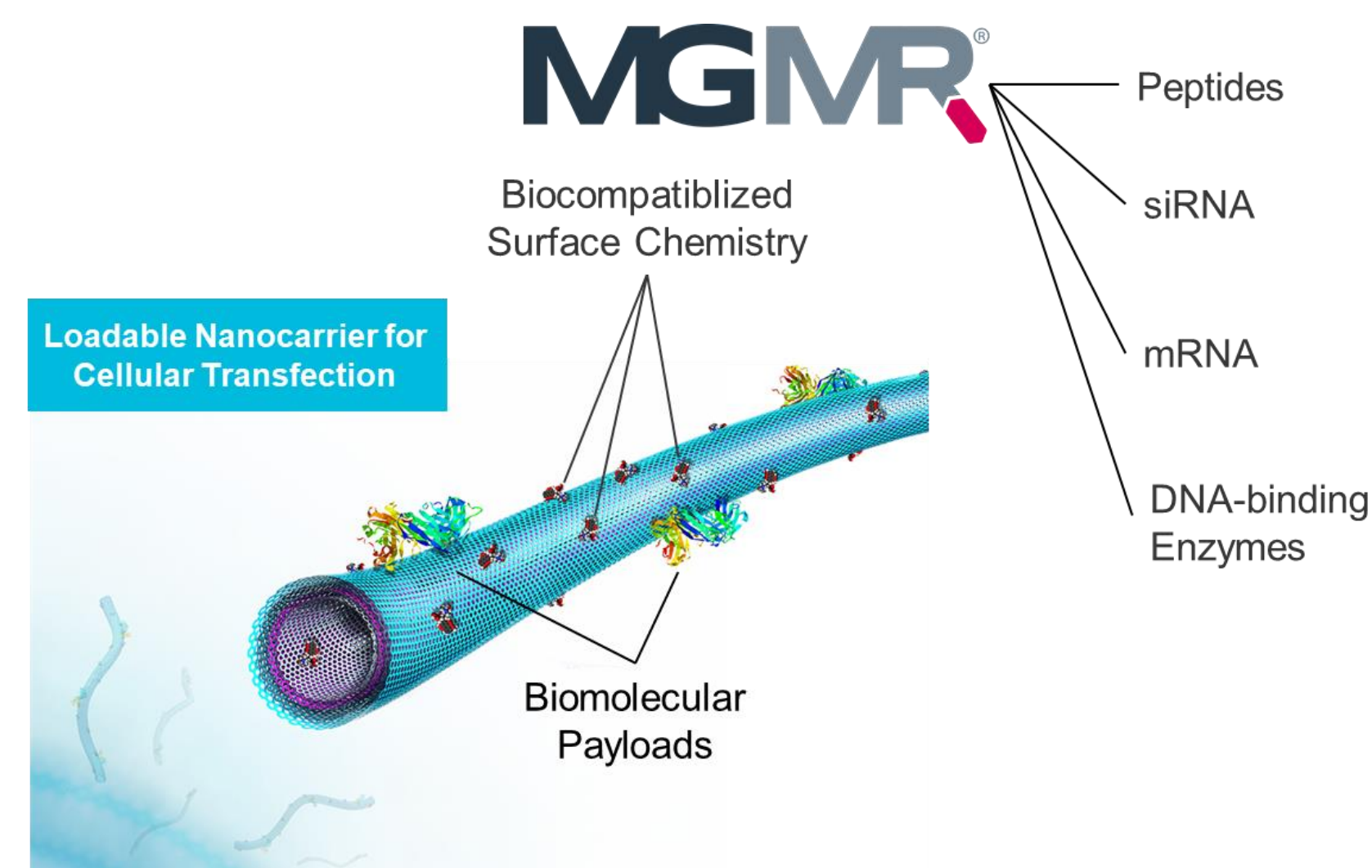


## Intracellular Transport Mechanism

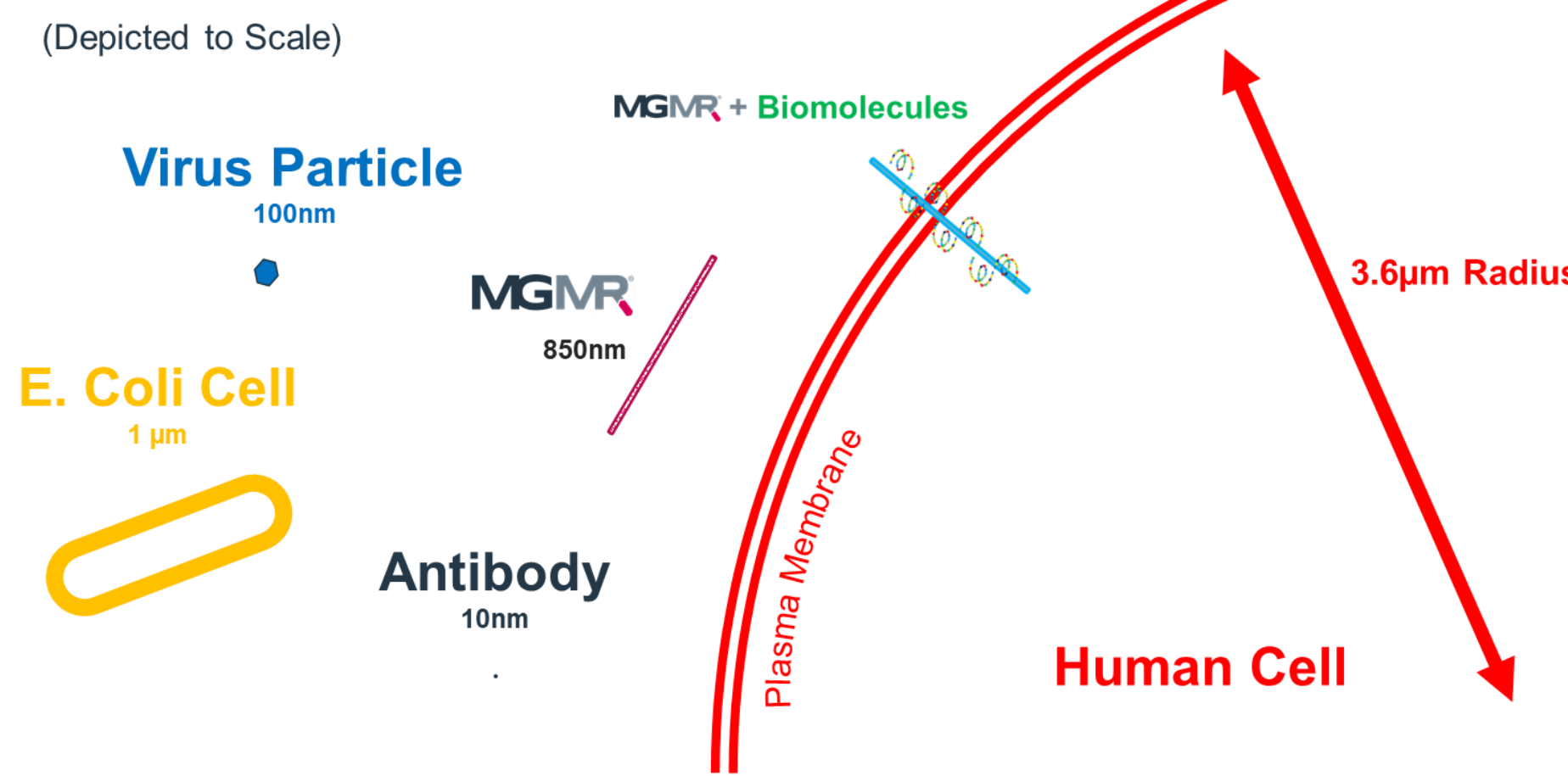

## MGMR<sup>®</sup> Supports T-cell Proliferation


### Dose-response Testing for Proliferation of Human T Lymphocyte-like cells (Jurkat)

- MGMR does not reduce cell proliferation relative to control
- (Growth Media, No Treatment)
- All groups indicated high viability
- Certain dosing-ranges of MGMR may improve cell proliferation and increase yield
- Protocol
  - Groups incubated with various concentrations of MGMR for 3 and 7 Days
  - Control = Standard Growth Media (RPMI 1640 + 15% FBS)
  - Cellular viability and quantification assessed by trypan exclusion assay
  - Stats: 95% CI

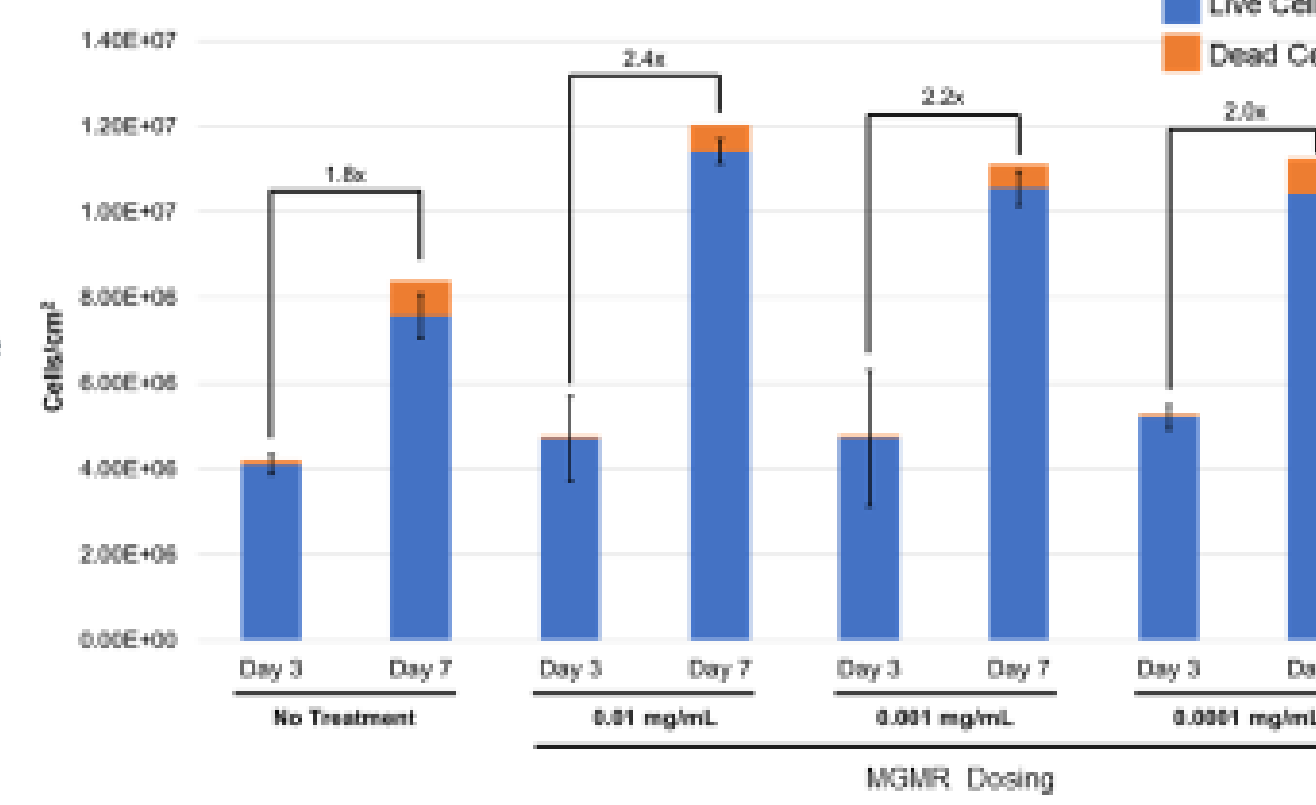

## Endocytotic Trafficking


- MGMR is transported into the cell by active transport
  - Unlike other transport systems, MGMR does not damage the cell's external casing (plasma membrane) in order to enter inside
- MGMR is initially encapsulated into larger compartments called vesicles - that traffic into the interior of the cell
  - MGMR is cargo agnostic - it is not prevented from getting inside cells by the size of its molecular cargo
- Once internalized, MGMR is released from vesicles into the intracellular environment
  - MGMR can deliver cargo to many different locations inside the cell

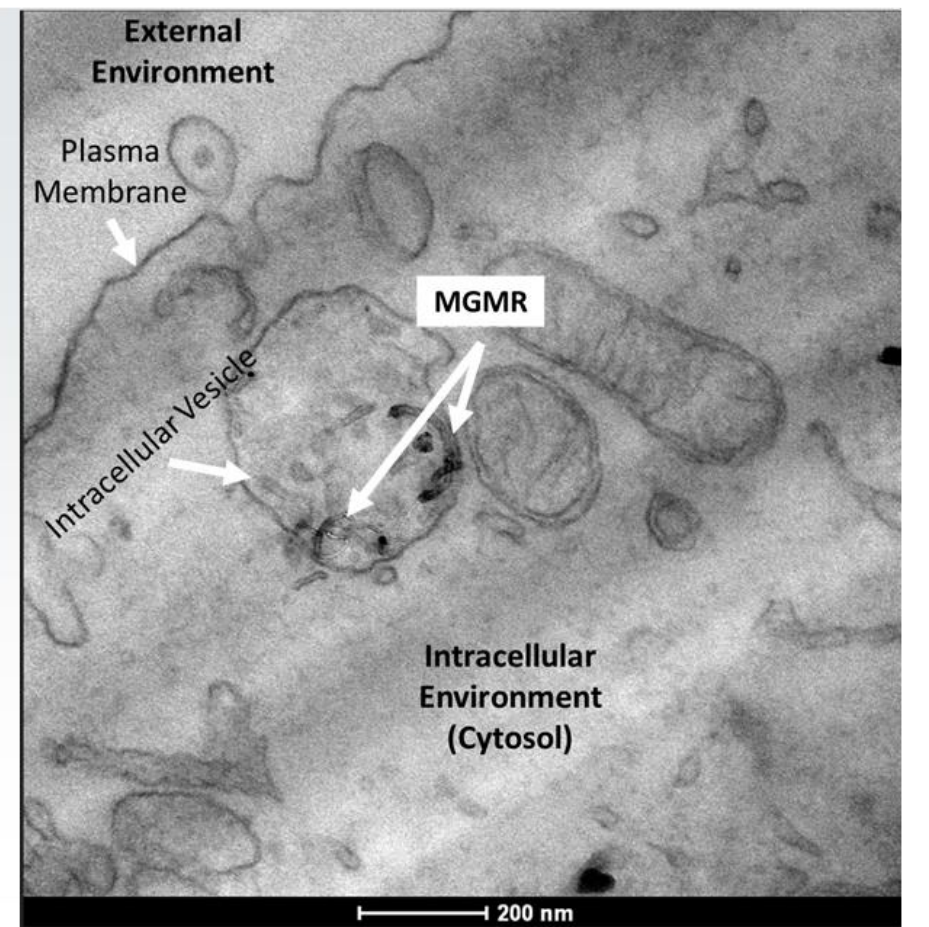



## MGMR<sup>®</sup> Peptide Delivery

MGMR demonstrated to enhance cell penetration of KLA—a cell “impermeable” 1.5kD protein<sup>1,2</sup>

- Proteins are powerful cell-regulators but very challenging for intracellular delivery
  - Do not efficiently cross the cell's membrane
  - Protein activity is very sensitive and easily diminished
- In this case study, MGMR binds to KLA with high affinity, transports it across the plasma membrane, and releases it—triggering apoptosis
  - MGMR+KLA peptide produced absolute apoptosis
  - MGMR alone has no effect (high tolerability)
  - KLA alone has no effect (benign externally)

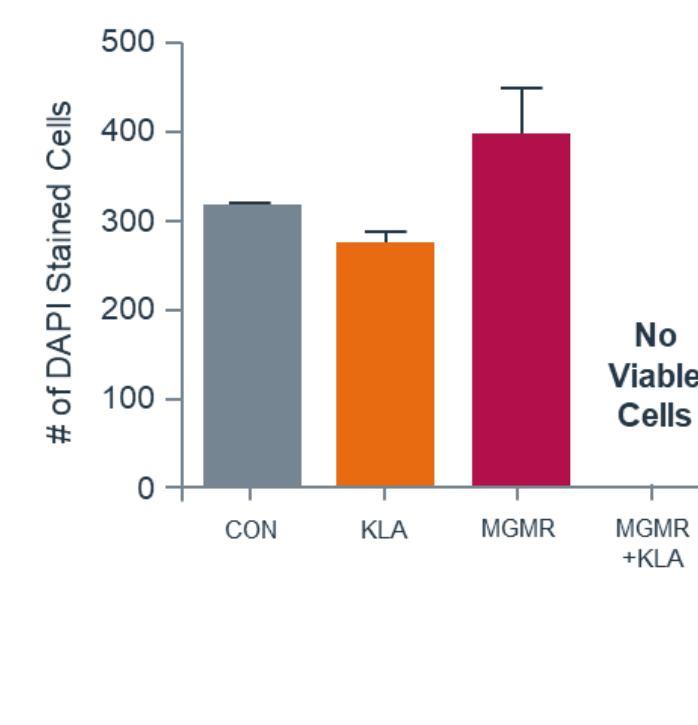


## MGMR<sup>®</sup> siRNA Delivery

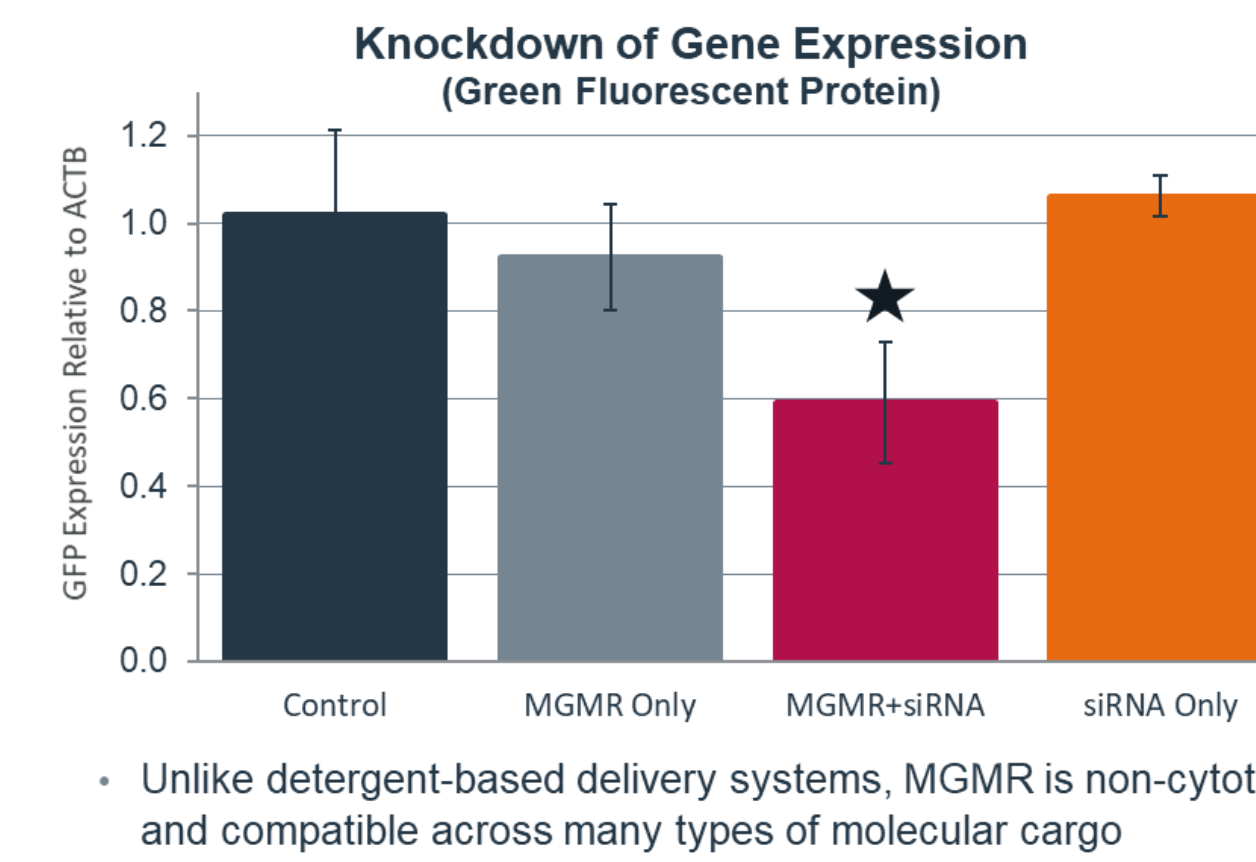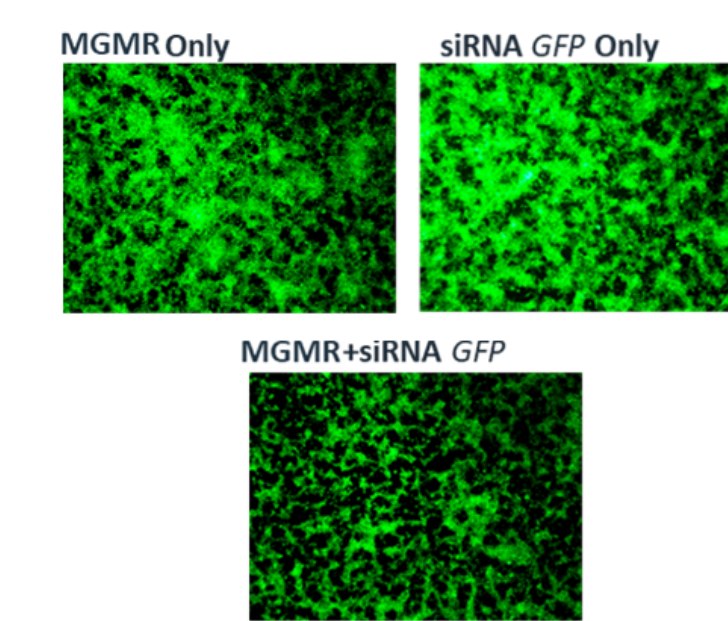


## MGMR<sup>®</sup> mRNA Delivery



### mRNA Delivery to T Lymphocytes (Jurkat)

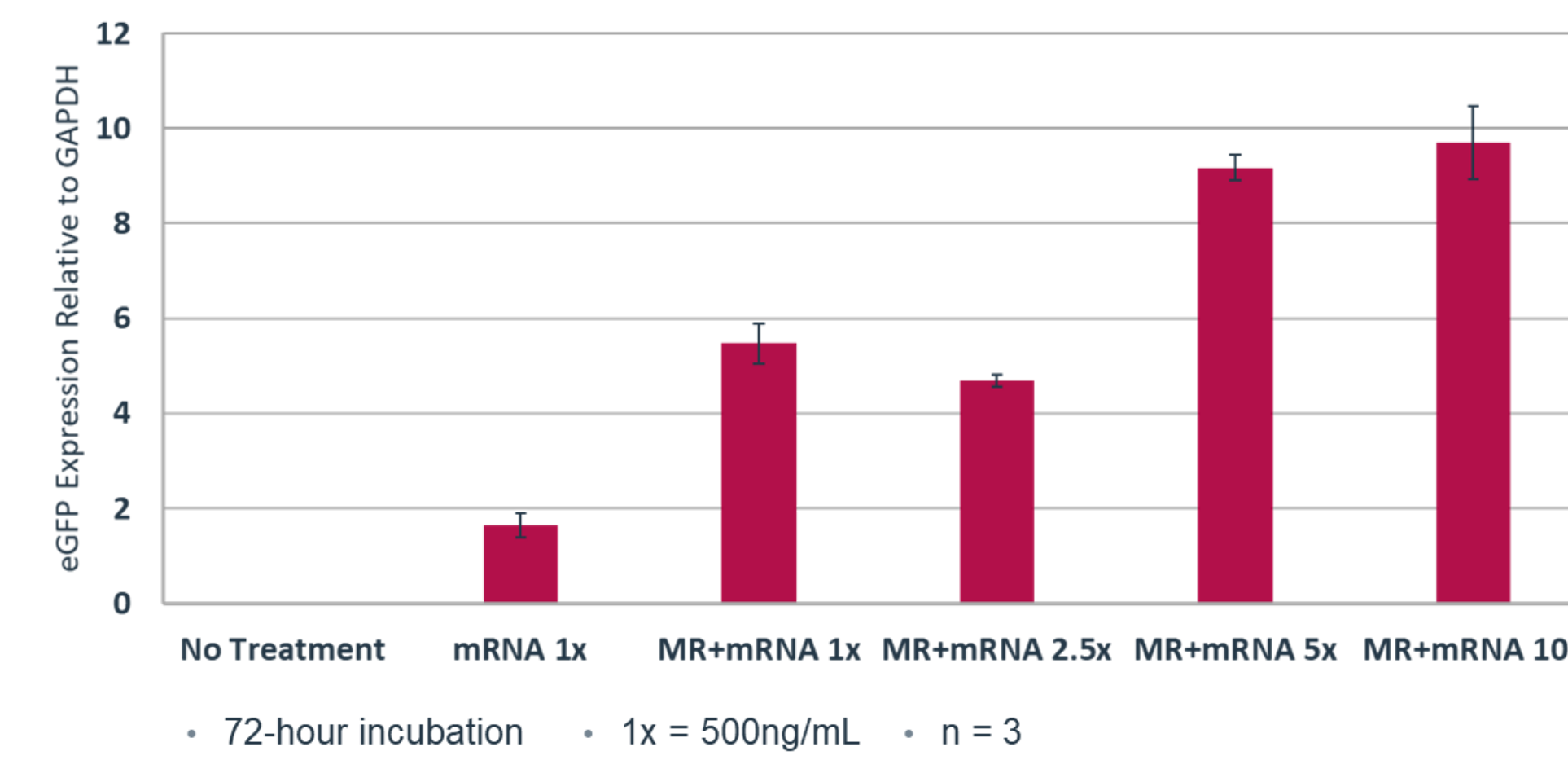


### mRNA Delivery to Monocyte-like Cells (U937)

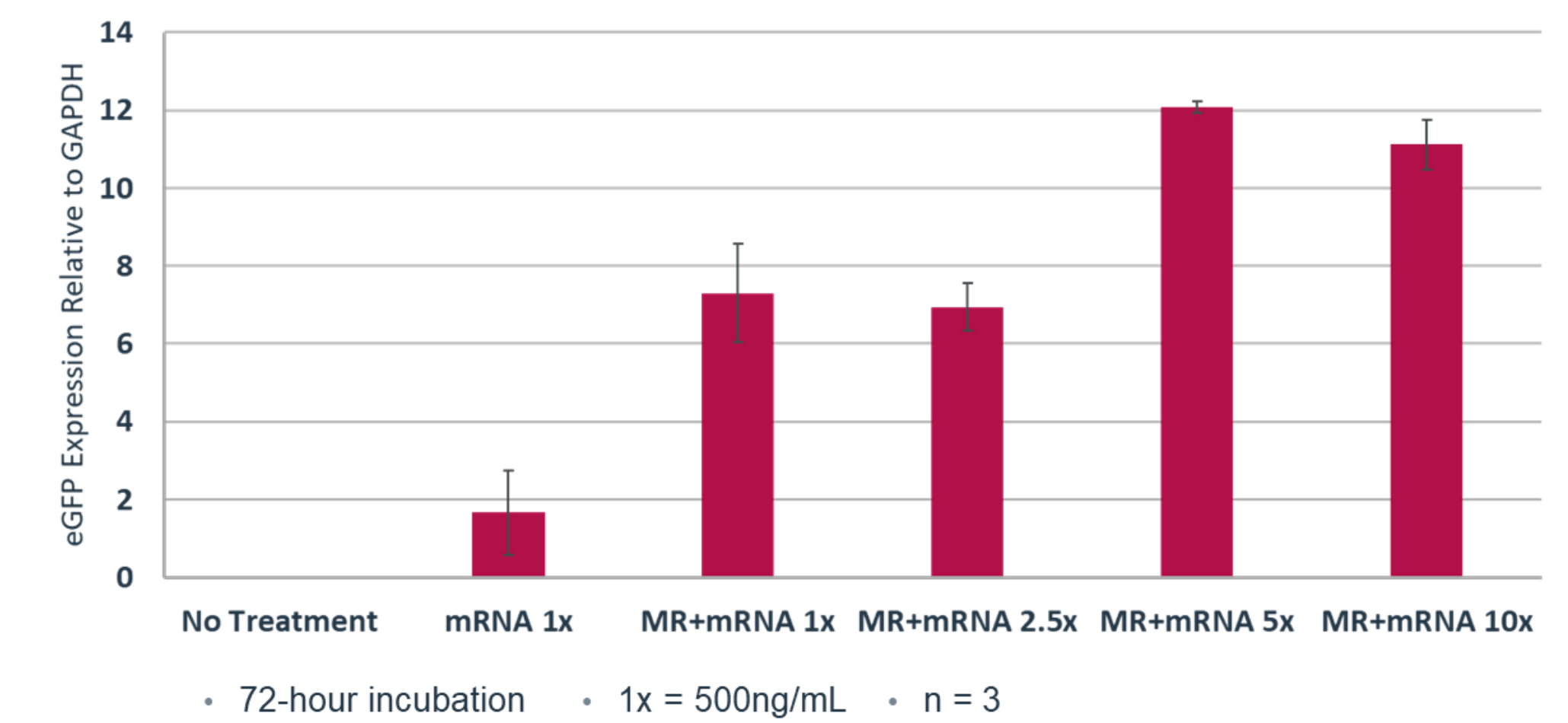


## MGMR<sup>®</sup> Nuclear Staining in Mesenchymal Stem Cells

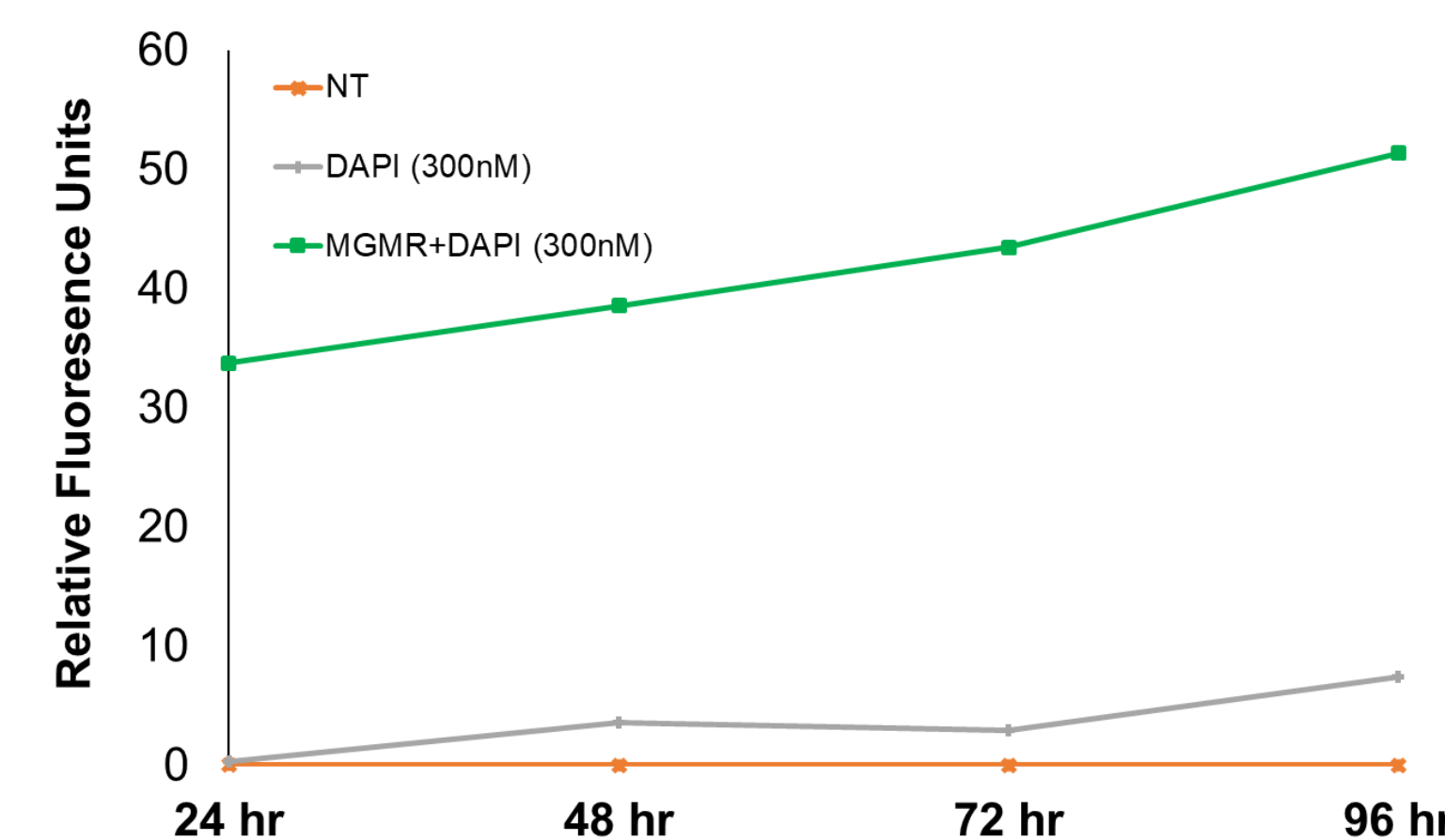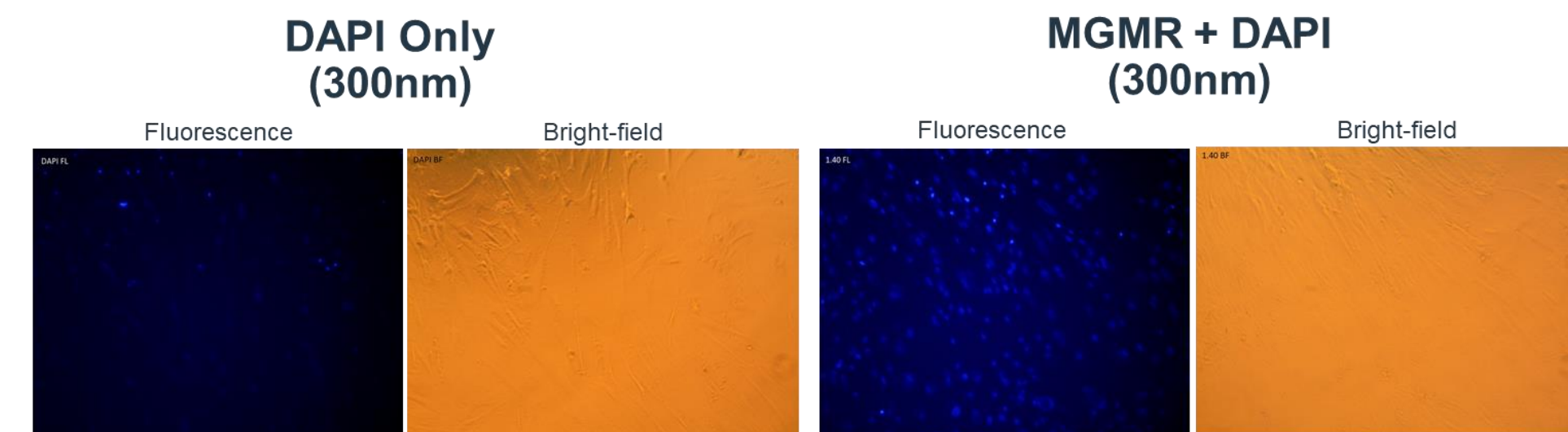

## MGMR<sup>®</sup> Delivery DNA-binding Enzymes in Stem Cells

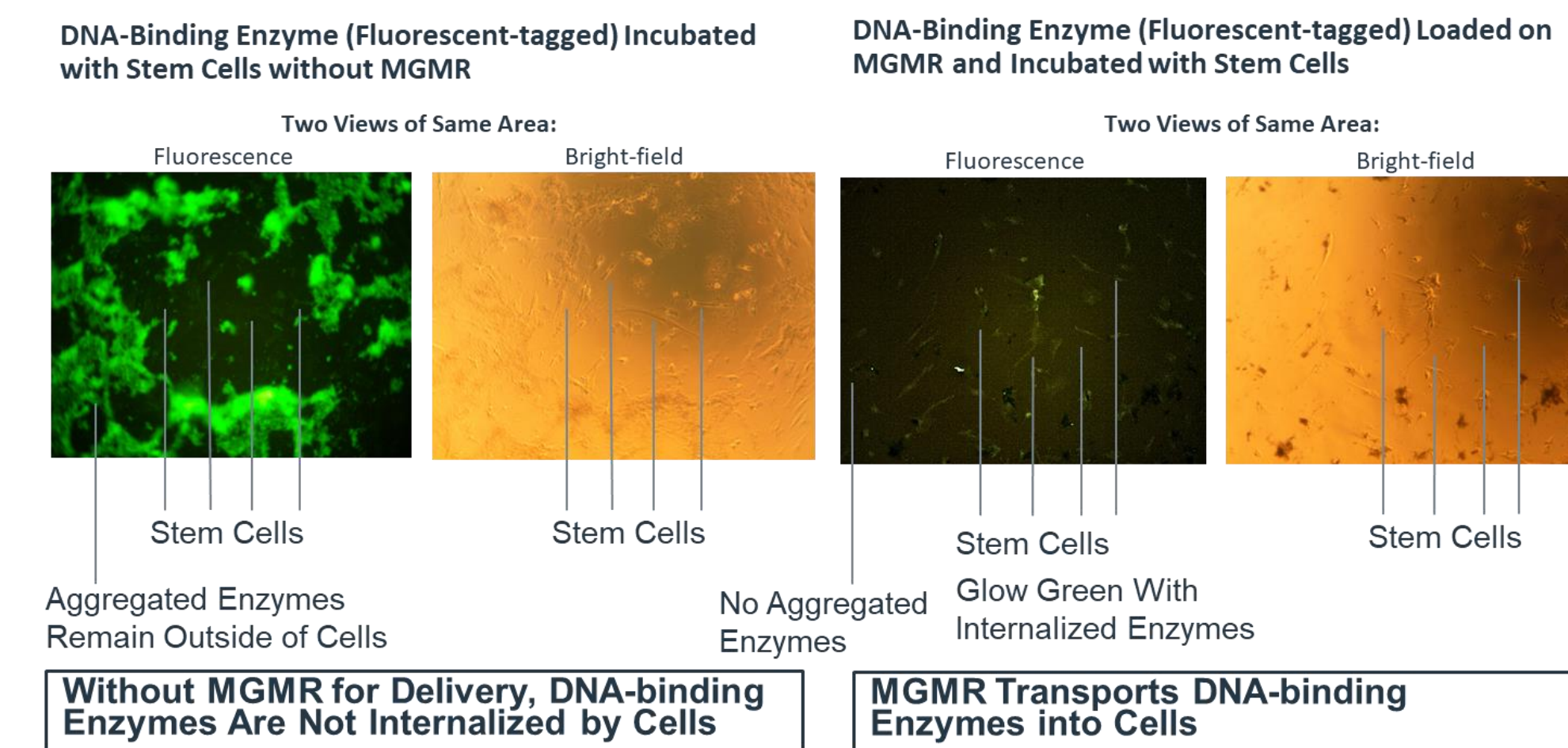

## MGMR<sup>®</sup> Delivery DNA-binding Enzymes in T-cells

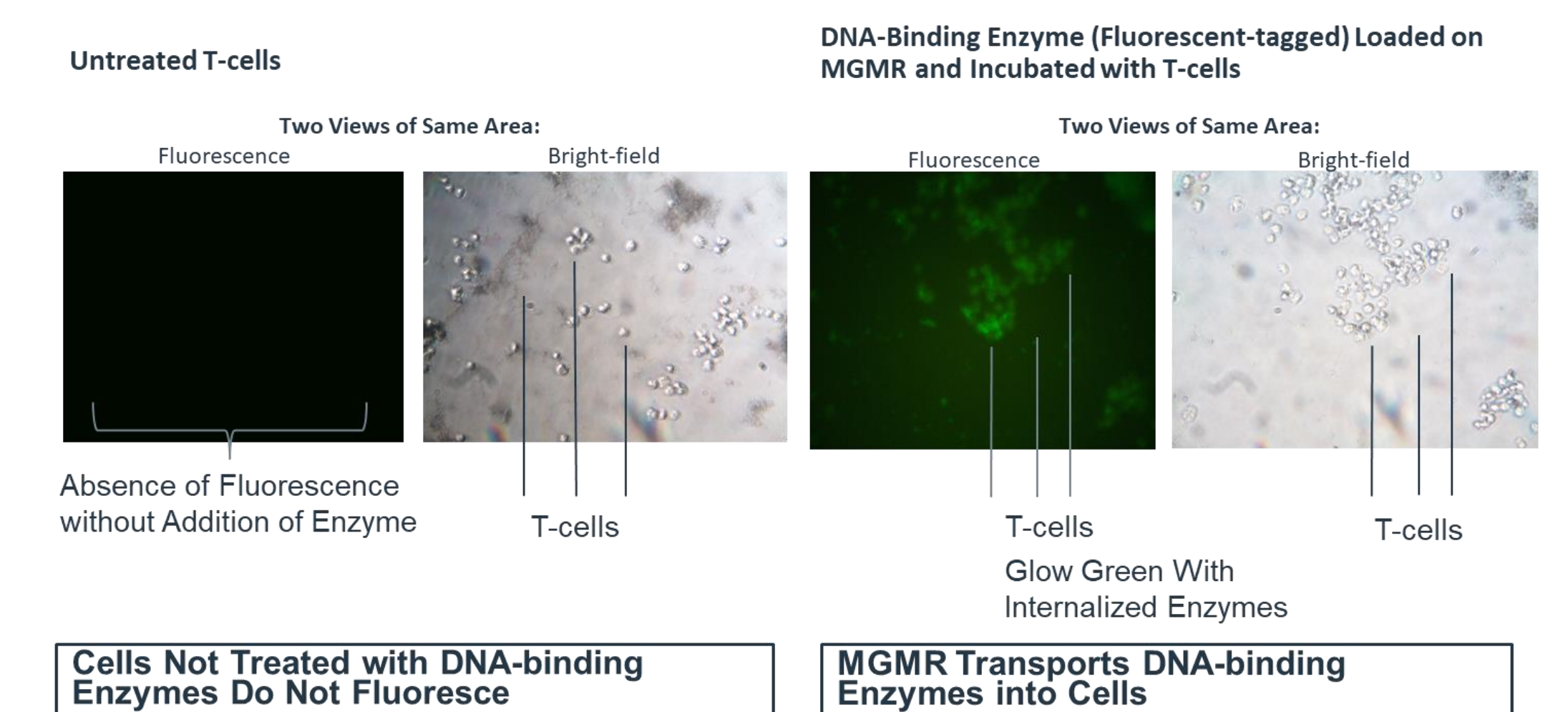

## MGMR<sup>®</sup> Next-generation Transfection Technology

Cargo	MGMR™	Viral Vectors	Polymers	Liposomes
Small Molecules	+	-	+	0
Peptides	++	+	+	0
Nucleic Acids	++	++	+	0
Proteins	++	+	+	0
Lipopolysaccharides	++	-	+	0
Log(P) Independent	+	+	-	-
Properties				
Non-immunogenic	+	-	+	+
Non-GMO	+	-	+	+
Simple Optimization	++	-	0	-
Scalable	++	+	0	-
Eukaryotes	+	-	0	0
Prokaryotes	+	-	0	-
Plants	+	-	0	-

**Key**  
++ Advantage  
+ Compatible  
0 Limited Compatibility  
- Non-Compatible

## How can MGMR<sup>®</sup> improve cell engineering?

- Eliminates** viral-mediated transfection
  - No risk of insertional mutagenesis
  - No viral toxicity and immunogenicity
  - Non-replicative
- Streamlines** manufacturing and quality processes
  - Non-biological transfection technology reduces variation
  - Readily extractable from final cell product
  - Eliminates need for new vector development – Single transfection vehicle for all cellular gene-editing systems
- Versatile** loading profile effective for intracellular delivery of:
  - Gene-editing complexes
  - Oligonucleotides
  - Peptides/Proteins
- Improves** cell yield
  - Reduced toxicity vs. viral or detergent-mediated transfection
  - Improved cell proliferation

## MGMR<sup>®</sup> Discover its Value in Your Process

Seamlessly Combine MGMR<sup>®</sup> with your Technology in 3 Phases

- Send Biopact Your Payload Molecule**
  - Provide Biopact with your gene-editing technology:
    - CRISPR
    - TALEN
    - ZFN
  - Biopact loads MGMR with your payload at **No Cost**
- Biopact Evaluates Performance**
  - Biopact will determine:
    - Loading efficiency on MGMR
    - Transfection efficiency in your target cell
  - Biopact reports performance of MGMR to you at **No Cost**
- Put MGMR<sup>®</sup> to the Test**
  - Order MGMR loaded with your molecule (\$10k-\$20k)
  - Readily sufficient for *in vitro*, *ex vivo* or *in vivo* testing
  - Evaluate MGMR loaded with your payload in your own studies

